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(URL <https://www.himac-science.jp>)

Keep this rotor instruction manual and the centrifuge manual in the file.

Angle Rotor for Refrigerated Centrifuge

R15A

Before using these rotors, please carefully read this instruction manual and the centrifuge instruction manual for its efficient operation and for your safety.

Keep this instruction manual for your reference and refer to it as required.

Contents

1. Specifications	_____	1
2. Tubes	_____	3
3. How to use	_____	4
4. Maintenance	_____	6
5. Packing list	_____	8
6. Spare parts	_____	8
7. Warranty of the rotor	_____	9
8. Decontamination	_____	10
9. Rotor retirement	_____	10
(Appendix) K Factors	_____	11

2021.03
S99864506

Original instruction manual

Copyright © 2021 Eppendorf Himac Technologies Co., Ltd.

All rights reserved. No part of this document may be reproduced or transmitted in any form or any means without permission from Eppendorf Himac Technologies Co., Ltd.

The names of actual companies and products mentioned herein may be the trademarks of their respective owners.

Safety Reminder

- Centrifuge rotors rotating at high speed have considerable potential for damage to personal properties if used improperly.
For safe and proper use of this rotor, carefully read the centrifuge instruction manual and this rotor instruction manual before use and observe the instructions.

 **WARNING :** and  **CAUTION :** notes are used to call your attention in this manual to prevent personal injury or damage to the rotor and the centrifuge.
These notes are defined as follows.

 **WARNING:** indicates a potentially hazardous situation which, if not avoided, could result in personal severe injury or possible death.
 **CAUTION:** indicates a hazardous situation which, if not avoided, could result in personal injury or severe damage to the instrument.

WARNING

- Never use any material capable of producing flammable or explosive vapors.
- Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.
- Never exceed the maximum speed of the rotor (mentioned on the rotor).
Always reduce rotor speed as instructed in this manual when rotor speed is limited due to sample density or kinds of tubes.
- Check the attached chemical resistance chart, and do not use any sample inapplicable to the rotor.
- If the centrifuge, rotor, or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the centrifuge, rotor, or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the centrifuge, rotor, or the accessory properly before requesting repairs from an authorized sales or service representative.
- It is your responsibility to sterilize and/or decontaminate the centrifuge, rotor, or parts properly before returning them to an authorized sales or service representative.

CAUTION

- Do not run the rotor without the rotor cover in position.
Tighten the cover-locking knob securely.
- Balance the tubes and the samples including the caps within the allowable imbalance of the rotor. Do not exceed the allowable imbalance.
- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Use the rotor tubes within their actual capacities.
- Do not use tubes that have exceeded their life expectancy. Failure to do so could result in damage of tubes and the rotor and the centrifuge. The life expectancy of tubes depends on factors such as the characteristics of samples, speed of the rotor used, and temperature.
Always check for deterioration and damage (cracks, deformation, and so on) on tubes before using them. Do not use the tubes if you find such a problem.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
Contact an authorized sales or service representative.

1. Specifications

1.1 Construction

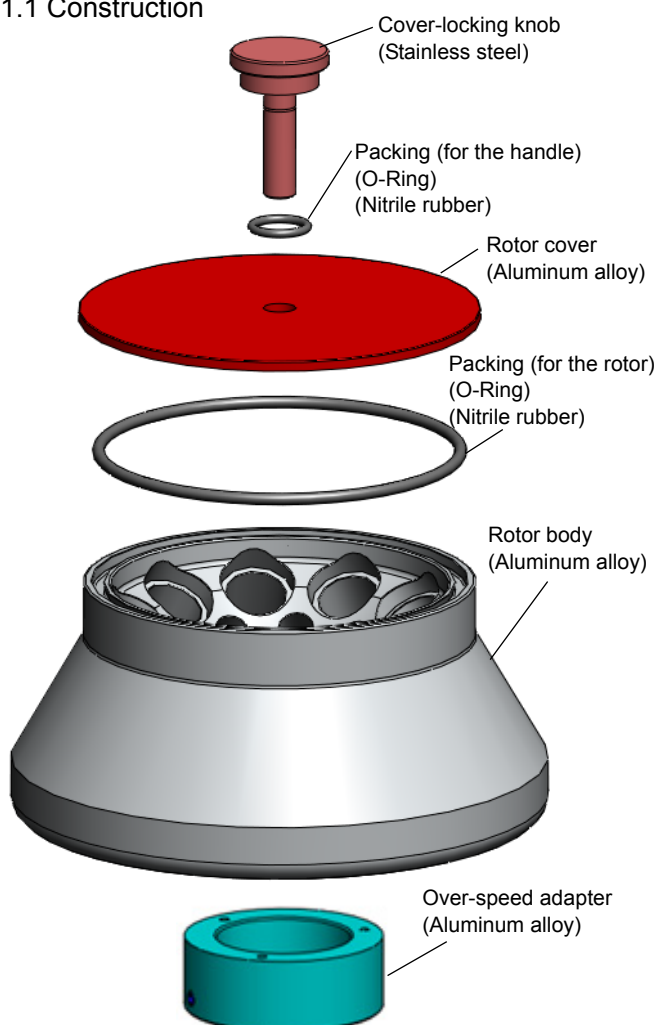
Figure1-1 Construction of the angle rotor

1.2 Specifications

Applicable centrifuge*1

(Some centrifuges are on sale in only Japan.
Visit our web site.)

Max. speed*2:	15,000 rpm
Max. RCF:	32,500 × g
Nominal rotor capacity:	50 mL × 10 tubes = 500 mL 15 mL × 10 tubes = 150 mL
Rotor Size:	φ 200 × 177.5 mm
Weight:	10.5 kg
Material:	Aluminum alloy
Rotor No.:	47

*1: For applicable centrifuges, refer to the “Applicable centrifuges (Rotors for high-speed refrigerated centrifuges (Part No. S998611))”.

*2: The maximum allowable speed varies with centrifuge models and so on. For further details, refer to the “Applicable centrifuges (Rotors for high-speed refrigerated centrifuges (Part No. S998611))”.

1.3 Cross-sectional view of rotor

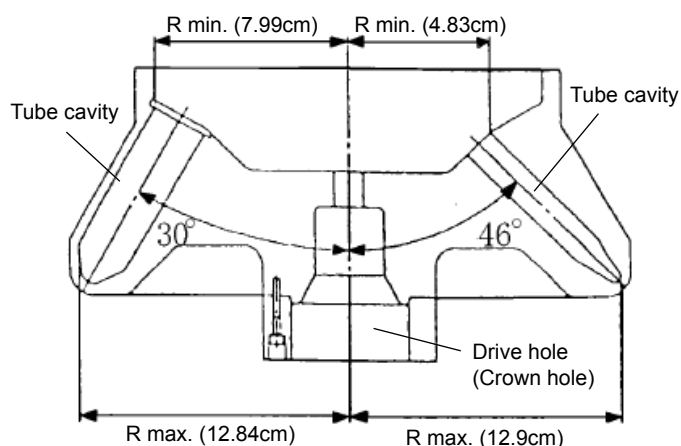
Figure1-2 Cross-sectional view of R15A rotor

1.4 Characteristic of rotor

Table1-1 Characteristics

(50mL Tubes)

Rotor speed (rpm)	RCF (×g)		K factor*
	Rmin. (7.99cm)	Rmax. (12.84cm)	
2,000	357	574	30,040
4,000	1,430	2,300	7,510
6,000	3,220	5,170	3,338
8,000	5,720	9,190	1,878
10,000	8,930	14,400	1,202
12,000	12,900	20,700	834
13,500	16,300	26,200	659
14,000	17,500	28,100	613
15,000	20,100	32,300	534

(15mL Tubes)

Rotor speed (rpm)	RCF (×g)		K factor*
	Rmin. (4.83cm)	Rmax. (12.9cm)	
2,000	216	577	62,210
4,000	864	2,310	15,552
6,000	1,940	5,190	6,912
8,000	3,460	9,230	3,888
10,000	5,400	14,400	2,488
11,000	6,530	17,500	2,056
12,000	7,780	20,800	1,728
14,000	10,600	28,300	1,270
15,000	12,200	32,500	1,106

*For details about K factor, refer to “(Appendix K Factors)”

Calculation formula

$$RCF = 1.118 \times 10^{-5} \times R \times N^2 (\times g)$$

R: Rotating radius (cm)
N: Speed(rpm)

1.5 Allowable speed of the rotor

WARNING :

Do not exceed the maximum speed of the rotor. The rotor speed should be limited depending on sample characteristic, kinds of tubes and centrifuge model.
Do not exceed the allowable speed of the rotor.

The maximum speed marked on the surface of the rotor cover is permitted under the conditions as follows.

-Average sample density : less than 1.2g/mL

-Standard tubes are used.

Depending on sample density, kinds of tubes and centrifuge model, the rotor speed may be limited as follows. Do not exceed the allowable speed of the rotor.

(1) Allowable speed for sample density

WARNING :

When sample density exceeds 1.2g/mL, calculate the allowable speed according to the following equation.

$$\text{Allowable speed (rpm)} = 15,000(\text{rpm}) \times \sqrt{\frac{1.2(\text{g/mL})}{\text{Average density of sample}(\text{g/mL})}}$$

(2) Allowable speed for combination of tube, cap, etc.

The rotor speed should be limited by combination of tube, cap, etc.

For details, refer to “2.1 Applicable tubes”.

Perform operation under the allowable speed, otherwise tubes may be broken during operation.

When using tubes on the market, perform operation under the allowable speed or the allowable RCF specified by the manufacturer. Otherwise tubes may be broken during the operation.

2. Tubes

2.1 Applicable tubes

CAUTION:

The tube cavities of this rotor are specifically designed for conical culture tubes.

Do not use round-bottom or self-standing culture tubes.

Use tubes which match the tube cavities of this rotor. Failure to do so, tube might be damaged and it might be hard to take out the tubes from the rotor.

- When using tubes on the market, perform operation under the allowable speed or the allowable RCF specified by the manufacturer and do not use any tubes whose allowable speed or RCF is not specified, otherwise the tubes may be broken during operation.
We will check the applicability of tubes upon request. Contact an authorized sales or service representative.
- Clean and sterilize tubes on the market according to the instructions of the manufacturer.
- Tubes on the market, himac 50TC tubes, and himac 15TC tubes should be discarded after one use.

Table2-1 Applicable tubes

Nominal capacity of the tube cavity of the rotor (mL)	Adapter		Tubes				Max. speed (Max. RCF)
	Name	Capacity × Qty	Part No.	Name	Actual capacity	Size (mm)	
50	—	—	S311187A	himac 50 TC tube(*1) (100 pcs./box)	42 mL	φ30 (*5) × 115	15,000rpm (32,300 × g)
	—	—	S307904A	himac 50 TC tube(*1) (300 pcs./box)	42 mL	φ30 (*5) × 115	15,000rpm (32,300 × g)
	—	—	(Marketed)	50mL PP culture tube (conical bottom)	42 mL	φ30 × 115	12,000rpm (20,670 × g)
15	—	—	S311184A	himac 15 TC tube(*1) (100 pcs./box)	15 mL	φ17 (*5) × 120	15,000rpm (32,500 × g)
	—	—	(Marketed)	15mL PP culture tube (conical bottom)	15 mL	φ17 × 120	11,000rpm (17,370 × g)
	Universal adapter (Eppendorf AG) (*2) (*3)	5 mL X 1	(Marketed)	Eppendorf Tubes® 5.0mL with screw cap (Eppendorf AG) (*4)	5.0mL	φ17 × 68	15,000rpm (22,400 × g)

PP: Polypropylene

(*1): Material of the tube: Polypropylene Material of the cap: Polyethylene

(*2): Order no. 5820 732.008 For the details, contact the maker of the adapter.

(*3): If it is hard to remove it from the tube cavity, use a tweezer etc. to pull it out of the tube cavity.

(*4): Order no. 0030 122.305 For the details, contact the maker of the tube.

(*5): Excluding the cap

The maximum allowable speed varies with centrifuge models and so on. For further details, refer to the “Applicable centrifuges (Rotors for high-speed refrigerated centrifuges (Part No. S998611))”.

The rotor speed is limited by the sample density and so on (see Section1.5).

Maximum speeds listed are guidelines only. Because of variances in user methodologies, no guarantee of performance is expressed or implied.

3. How to use

3.1 Preparation of tubes

Check that the tubes are free from crack or deformation. Do not use remarkably deformed or cracked tubes.

WARNING :

- This rotor and the centrifuge are not explosion-proof. Never use explosive or flammable samples.
- For safety, there are limits on the use of bio-samples which require bio-isolation, such as pathogenic germs and DNA recombination, as well as RI substance in centrifuges. Perform strict safety controls when separating samples containing these substances.

CAUTION:

- Do not use any sample inapplicable to the tubes and the caps in use referring to the attached "Chemical resistance chart". Otherwise, the tubes and the caps may be degraded.
- Chemical resistance and the strength of tubes varies with speed, temperature, and so on. Before using sample, fill tubes with water, buffer solution, etc. instead of sample and run them at the intended speed, temperature, and so on to ensure that there is no abnormality.
- Do not exceed the specified sample capacity of the tubes, otherwise sample leakage or deformation of tubes may occur.
- Balance the sample within the allowable imbalance including the weight of the tubes, caps, and adapters. Excessive imbalance can cause damage to the centrifuge and the rotor.
- Use the samples that are same in compositions and same in density when placing them symmetrically. Otherwise the precipitation levels may be different by centrifugation and such operation may increase the level of imbalance, and it may result in damage to the centrifuge and the rotor.
- Do not exceed the actual capacity specified in the manufacturer. Otherwise the sample leakage may occur during operation.
- Use the tubes that are same products. Otherwise, the tubes may be damaged due to their different form or the imbalance operation may occur, and it may result in the damage to the centrifuge and the rotor.
- Use the tubes at a temperature between 4°C and 25°C. Otherwise, tubes may be broken or deformed during operation.
- Before using tubes on the market, test them by filling with water instead of sample, and run them at the intended speed to ensure that there is no abnormality.

Inject the samples into the tubes and cap the tubes properly. Do not exceed the specified sample capacity of the tubes, otherwise sample leakage or deformation of tubes may occur.

Balance the tubes that are symmetrically arranged. The allowable imbalance of this rotor is within 2 g.

3.2 Setting of tubes into rotor


CAUTION:

Do not load only one tube nor arrange the tubes asymmetrically with respect to the drive shaft, otherwise the centrifuge or the rotor may be damaged due to excessive imbalance operation.

Check that the tube cavities of the rotor are free from foreign substances. Set the balanced tubes symmetrically with respect to the drive shaft.

Pre-cool the rotor when the operating temperature of the rotor is lower than the room temperature or pre-heat the rotor when the operating temperature of the rotor is higher than the room temperature before setting the tubes in the rotor.

Examples of incorrect tube setting are shown below.

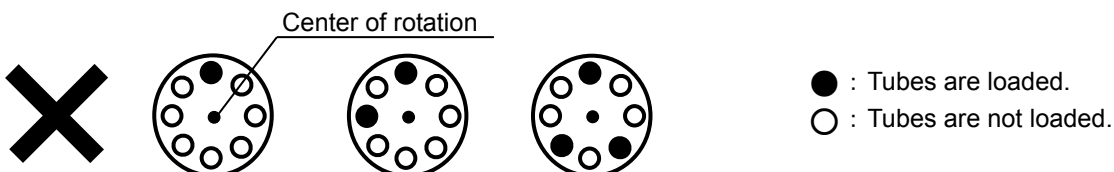3.3 Mounting the rotor cover


CAUTION:

- Be sure to mount the rotor cover to the rotor body and tighten the cover locking knob securely before operation, otherwise the rotor or the rotor cover may be removed during operation and result in damage to the centrifuge and the rotor.
- Do not mount the rotor cover into the rotor while the rotor is placed in the rotor chamber. Doing so might cause the rotor cover to be tightened insufficiently.
Do not remove the rotor cover from the rotor while the rotor is placed in the rotor chamber. Doing so might cause the parts of the rotor cover to be damaged.

- (1) Apply silicone grease (vacuum grease) (standard accessory of the centrifuge) to the packing (for the rotor) and then fit the packing (for the rotor) in the groove of the rotor body. Replace the packing (for the rotor) with new one if damaged or deteriorated.

If you neglect to apply silicone grease (vacuum grease), it might be hard to mount or remove the rotor cover. When it is hard to mount or remove the rotor cover, apply silicone grease (vacuum grease) to the packing (for the rotor).

- (2) Apply lubricant for screw (Part No. 84810601) to the threaded portion of the cover locking knob. Put the rotor cover on the rotor body. Turn the cover locking knob to screw in the rotor body securely.

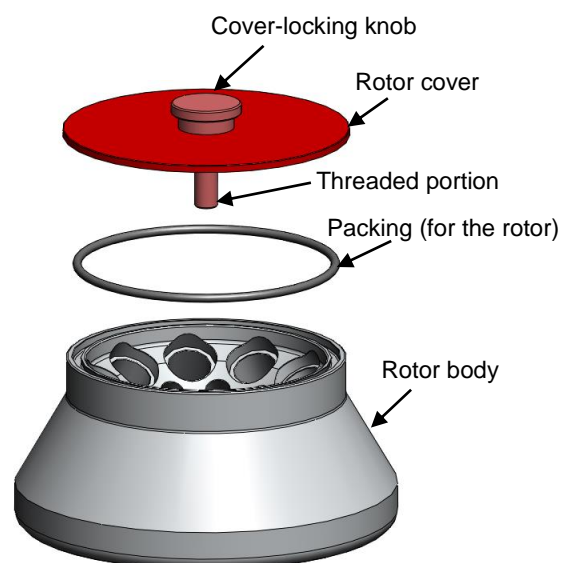3.4 Operation

Mount the rotor onto the drive shaft of the centrifuge gently and securely. Check that the rotor cover is mounted and the cover-locking knob is securely tightened. Perform operation according to the instructions of the centrifuge instruction manual.

3.5 Taking out samples

After the operation, gently take out the rotor from the centrifuge and remove the rotor cover by turning the cover locking knob. Take out the tubes from the rotor and then take out the sample in a proper manner.

4. Maintenance

4.1 Corrosion resistance of rotor

WARNING:

Check the attached chemical resistance chart, and do not use any sample inapplicable to the rotor.

This rotor is made of aluminum alloy. The rotors made of aluminum alloy have high corrosion resistance and they are covered with an anodic oxidation coating. However, use of inapplicable chemicals can corrode these rotors and decrease the strength. Use samples that will not affect the rotor referring to the attached chemical resistance chart.

4.2 Maintenance of rotor

After the operation, properly maintain the rotor to prevent corrosion that can cause the rotor breakage.

- Normal maintenance

CAUTION:

Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

Wash the rotor with tap water or a dilute solution of neutral detergent and rinse it out with distilled water. Wipe the rotor with a soft cloth. Turn the rotor upside down with the rotor cover removed to dry it well. Check that the rotor is completely dried and then put a light coat of silicone grease (vacuum grease) (standard accessory of the centrifuge) on the rotor. Store the rotor in a dry place.

- Maintenance after use of a corrosive sample

Wash the rotor with tap water immediately after the operation. Then perform the normal maintenance.

- Maintenance when foreign substances are adhered to the rotor

WARNING :

Do not allow the rotor temperature to rise over 100°C. Otherwise, the rotor can be brittle.

CAUTION:

Use a neutral detergent having a pH between 5 and 9, otherwise the rotor can be discolored or corroded.

Soak the rotor in warm water for one or two hours and wash the inside of the tube cavities and the crown hole at the rotor bottom with a soft brush to remove foreign substances. Do not forget to remove foreign substances adhered in the crown hole at the rotor bottom, otherwise the rotor and the drive shaft may be damaged.

4.3 Sterilizing rotor

- Sterilize the aluminum alloy rotors according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin).

WARNING :

- Never sterilize the rotor by autoclaving or boiling. Otherwise, the rotor will be brittle. It is very dangerous.

CAUTION:

- Do not dip the rotor in the formalin (3%) solution more than 2 hours.

4.4 Inspecting rotor


Periodically check that the rotor is free from corrosion every 100 hours use. Especially be careful of the tube cavities of the rotor and the crown hole at the rotor bottom because the rotor can be brittle if these portions are corroded. If the rotor surface is discolored, dented or cracked, the rotor is corroded. Do not use such a corroded rotor and contact an authorized sales or service representative immediately for further inspection.

5. Packing list

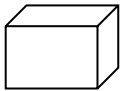
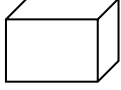
Check the parts and accessories of the rotor when the rotor is delivered. Contact an authorized sales or service representative if there are a missing item and inferior goods.

No.	Name	Qty	Remark
①	Rotor	1	—
②	Packing (for the handle) (O-Ring)	1*	—
③	Packing (for the rotor) (O-Ring)	1*	—
④	himac 50TC tube	1	50TC Tube: 100 pcs. 50TC Cap: 100 pcs.
⑤	himac 15TC tube	1	15TC Tube: 100 pcs. 15TC Cap: 100 pcs.
⑥	Rotor instruction manual	1	Part No. S998645

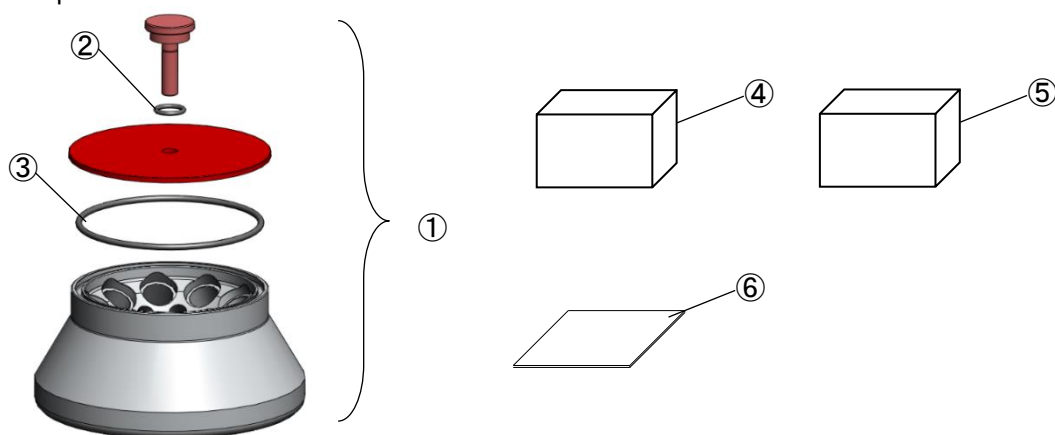*: the quantities of the parts that are mounted to the rotor

6. Spare parts

Figure	Part No.	Name	Qty
	84521016	Packing (for the handle)(O-Ring)	5
	84521047	Packing (for the rotor)(O-Ring)	3
	S311187A	himac 50TC tube	100
	S311184A	himac 15TC tube	100
	84810601	Lubricant for screw	1
	483719	Silicone grease(vacuum grease)	1
	S998645	Rotor instruction manual	1

7. Warranty of the rotor

●Warranty for Refrigerated Centrifuge Rotors

Eppendorf Himaс Technologies warrants its refrigerated centrifuge rotors under the terms and conditions set forth below.

1. Scope of Application

This warranty applies to:

- (1) failure or damage to the rotor due to rotor defects, in materials or workmanship, that are attributable to Eppendorf Himaс Technologies, and
- (2) failure or damage to an Eppendorf Himaс Technologies refrigerated centrifuge due to defects in a rotor covered by the warranty.

Note: This warranty covers the rotor or refrigerated centrifuge only, and Eppendorf Himaс Technologies shall not be liable for damage to samples or other items that may arise from failure of or damage to the rotor or refrigerated centrifuge.

2. Warranty Period

This warranty is valid for seven years from the date of delivery of the rotor by Eppendorf Himaс Technologies or an Eppendorf Himaс Technologies dealer.

3. Notes

- (1) This warranty extends only to the original buyer from Eppendorf Himaс Technologies or the original buyer from an Eppendorf Himaс Technologies dealer. Rotors whose ownership passes from the original buyer to a third person due to resale, transfer or the like are not covered by this warranty. Neither does this warranty cover cases where the ownership of a rotor is reacquired subsequently by the original buyer.
- (2) The warranty for the rotor proper (warranty period and terms) shall continue to apply in cases where rotor parts and the like are additionally purchased.
In such cases, the warranty period shall continue to be measured from the date of delivery of the rotor. If such parts are subject to a restricted warranty period, this warranty shall only cover failure or damage involving such parts that occurs both within the rotor warranty period and within the warranty period for the parts concerned.
- (3) This warranty does not cover any of the following, even during the warranty period:
 - ① Failure or damage due to abuse or misuse.
 - ② Failure or damage due to operation or maintenance in a manner contrary to the instructions in the operator's manuals for the rotor and refrigerated centrifuge.
 - ③ Failure or damage due to disassembly or modification without the permission of Eppendorf Himaс Technologies.
 - ④ Failure or damage due to use of a bucket, adapter, tube, bottle, or the like, other than those approved by Eppendorf Himaс Technologies for use with the product.
 - ⑤ Failure or damage due to disaster such as fire or earthquake.
 - ⑥ Failure or damage due to use with a refrigerated centrifuge not of Eppendorf Himaс Technologies manufacture without the permission of Eppendorf Himaс Technologies.
 - ⑦ Failure or damage due to use with a refrigerated centrifuge that has been modified without the permission of Eppendorf Himaс Technologies.
 - ⑧ Failure or damage due to a failed or damaged part (including a tube, bottle, adapter, etc) not submitted to Eppendorf Himaс Technologies.
 - ⑨ Replacement of consumable items.
 - ⑩ Parts subject to a restricted warranty period that has expired.

●Other Matters

Rotors that are delivered to Eppendorf Himaс Technologies for inspection or repair must be sterilized and cleaned beforehand.

Be aware that Eppendorf Himaс Technologies will sometimes decline to carry out inspection or repair if a rotor or other related equipment is or could be contaminated as a result of use without protection or under inadequate safety management in a radioisotope facility or a laboratory of the level P2 or higher.

8. Decontamination

WARNING:

- If the centrifuge, rotor or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the rotor or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the rotor or the accessory properly before requesting repairs from an authorized sales or service representative. Note that we cannot repair the centrifuge, rotor or the accessory unless sterilization or decontamination is completed.
- It is your responsibility to sterilize and/or decontaminate the rotor or parts properly before returning them to an authorized sales or service representative. In such cases, copy the attached decontamination sheet and fill out the copied sheet, then attach it to the item to be returned. We may ask you about the treatment for the rotor or parts if the decontamination is checked and judged as insufficient by us. It is your responsibility to bear the cost of sterilization or decontamination. Note that we cannot repair or inspect the rotor or the accessory unless sterilization or decontamination is completed.

9. Rotor retirement

After many years of use, there will be inevitably some corrosion or stress corrosion. At some points, the combination of such damage and metal fatigue could make the rotor vulnerable to a failure.

Although a rotor may appear to be in a good condition, you should follow the rotor retirement recommendation shown below.

Rotor	Material	Retire After Years
Refrigerated centrifuge rotors	Aluminum alloy	15
	Stainless steel	

(Appendix) K Factors

- Calculation of separation time

The K factor can conveniently be used which makes it possible to estimate the settling time directly from Svedberg unit(S).

Namely, the time taken for particles of 1S to settle from Rmin. to Rmax. is K factor.

The K factor can be obtained from the formula show below.

$$K = \frac{\ln(R \max.) - \ln(R \min.)}{\omega^2} \cdot \frac{10^{13}}{3600}$$

ω : Angular velocity $\omega = 2\pi/60 \times N$

N : Speed (rpm)

Rmin. : Distance from axis of rotation to top of solution (cm)

Rmax. : Distance from axis of rotation to bottom of tube (cm)

Using the K factor, the settling time t (in hours) is expressed as follows:

$$t = \frac{K}{S}$$

For example, the K factor at the maximum speed of 15,000rpm of Model R15A rotor is 534. The time taken for particles of 100 S to be settled from Rmin. to Rmax. using this rotor is as follows:

$$t = \frac{534}{100} = 5.3 \text{ hours}$$

The time taken for settling of particles of 100 S by the same rotor at a speed of 15,000 rpm is:

$$K = \left(\frac{15,000}{12,000} \right)^2 \times 534 = 834 \quad t = \frac{834}{100} = 8.3 \text{ hours}$$

Eppendorf Himac Technologies Co., Ltd.

1060, Takeda, Hitachinaka City
Ibaraki Pref., 312-8502 Japan

URL: <https://www.himac-science.com>

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Copyright © 2021 Eppendorf Himac Technologies Co., Ltd. All rights reserved.